
MNova Installation and Activation Guide

The Chemistry Institutes of the LUH hold a Campuse license of the software MNova for processing of NMR data which includes the software packages qNMR and Reaction Monitoring. The campus license files are updated once a year to the latest version.

The license files have to be activated and verified by a connection to the license server which can be accessed from the university network. If no LAN connection to the university network exists, a VPN connection is required. The WiFi connection via eduroam is not sufficient for the license activation. For institutes other than the OCI a VPN connection may also be required.

Instructions for setting up a VPN network are given by the [LUIS](#). The activated license is valid for 90 days during which MNova can be used offline. The license file has to be verified again by connecting to the license server at the latest each 90 days.

You can download the latest MNova version from their [website](#). Please make sure that you are in the LUH and/or VPN network upon starting MNova the first time.

You can download the license files from the homepage of the NMR department or from this [link](#).

The license files can be installed in two different ways:

- (1) Unzip the license files and copy them into the directory (C:/Program Files/)Mestrelab Research S.L/MestReNova/licenses. The license file will be activated automatically upon the first start.
- (2) Unzip the license files into any folder on your hard disk. At your first start of MNova you will be asked for the license files. Click install and browse for a license file. This step has to be repeated for every license file.

Troubleshooting

1. Check your VPN connection.
2. Download the latest license files.
3. Check the license manager in MNova for the installed license files. You find the icon in the upper right corner besides the "Options" menu.
3. If this does not solve your problem please contact the NMR staff.